

Encapsulator RX Directions

FOR WATER-BASED SPILLS QUICKLY ABSORBS: ACIDS AND CAUSTICS

Directions:

Use Encapsulator Rx to form a containment dike around the spill. Then cover the spill completely. Add more as needed. The absorbed spill will range from a soft to a hard gel. The absorbed spill may then be picked up and removed from the spill site.

Encapsulator Rx will absorb liquids and gasses.

Encapsulator **Rx** will not promote Off-Gassing.

Depending on the size & make-up of the spill: Neutralize and dispose of accordance with federal, state and local regulations.

CAUTION: Do not spray with water, it may cause slipping. If dusting should occur, do not breathe the dust for extended periods of ime. With any liquid corrosive spill the use of gloves, goggles, and respirator is recommended.

Keep lid tightly closed when not in use.